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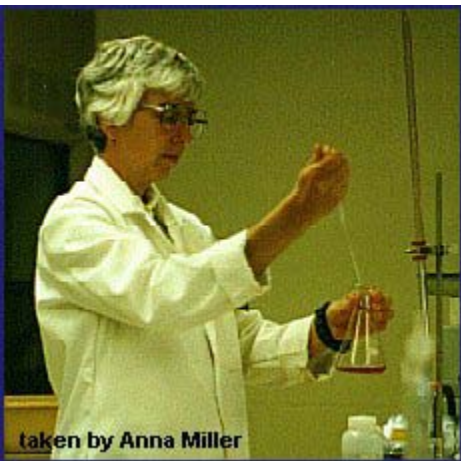
Hood sash  
set above  
18 inches.

- Excess storage of chemicals.
- Exhaust slots blocked.
- Containers stored within six inches of face of hood.

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- **Chemical Fume Hoods**
- Responsibility: EHS.
- Testing shall occur with HVAC systems operating at 100%. Qualified and trained personnel shall conduct testing. At a minimum, the following test procedures will be followed:
- Visual inspection for hood damage, modifications or congestion.
- Face velocity measurements at the designated working sash position. (To pass, the average velocity must be plus or minus 10% of the design rating at the standard operating sash opening. For most chemical fumehoods at the University, standard operating sash opening is 18 inches sash height.)
- Chemical fumehood alarm test.
- University of North Carolina at Chapel Hill
- 02 June 2011 0-00-7
- Tested chemical fumehoods will have EHS labels affixed stating the following: test date; initials of inspector, and face velocity readings at the standard operating sash opening as indicated on the test label.
- For chemical fume hoods found to be above or below 10% of the design rating below at the standard operating sash opening, Facilities Services will be contacted for airflow adjustments. (this does not apply to low flow chemical fume hoods) A low flow hood is one that has had the exhaust volume reduced by operating through a smaller sash opening. This type of hood does not require the containment be the same with the sash full open for setup as it is for usage. Energy savings can parallel that of low velocity hoods, but proper sash position for usage or setup must be managed. EHS will post hoods: "Out of Service/Do Not Use." Hoods that are posted "Out of Service/Do Not Use" may need to be emptied and cleaned as determined by the EHS Safety Clearance procedure to prevent chemical exposures for lab occupants and personnel who will repair the hood system.